## BG-Power600k Power Supply quick start manual

### Quick start

#### Setting voltage, current or power limits

The display will now flash for the set voltage limit. Using the SCROLL KEY to setting voltage limit and then switch to suitable digit using the SWITCH KEYS. Using the SCROLL KEYS again to select the desire values for the run. Confirm with ENTER KEY. Repeat above procedures for limiting current and power.

#### Setting time limit

The default time is 999 minutes. In case changing the time limit, switch to suitable digit using the SWITCH KEYS and select the desire values with the SCROLL KEYS. Press RUN and PAUSE KEY to run.


### Program

#### Saving a program

Choose a blank program and set the voltage, current and power limit as described above. Press RUN and PAUSE KEY will run the parameters. By the end, the parameters will be saved to the blank program. The instrument can save up to twelve programs.

#### Choosing a program

Using the SCROLL KEYS to select a suitable program and press ENTER to retrieve the parameters. Press RUN and PAUSE KEY to run. The parameters can also be changed using the procedures described above, and the changes will be saved to the program.

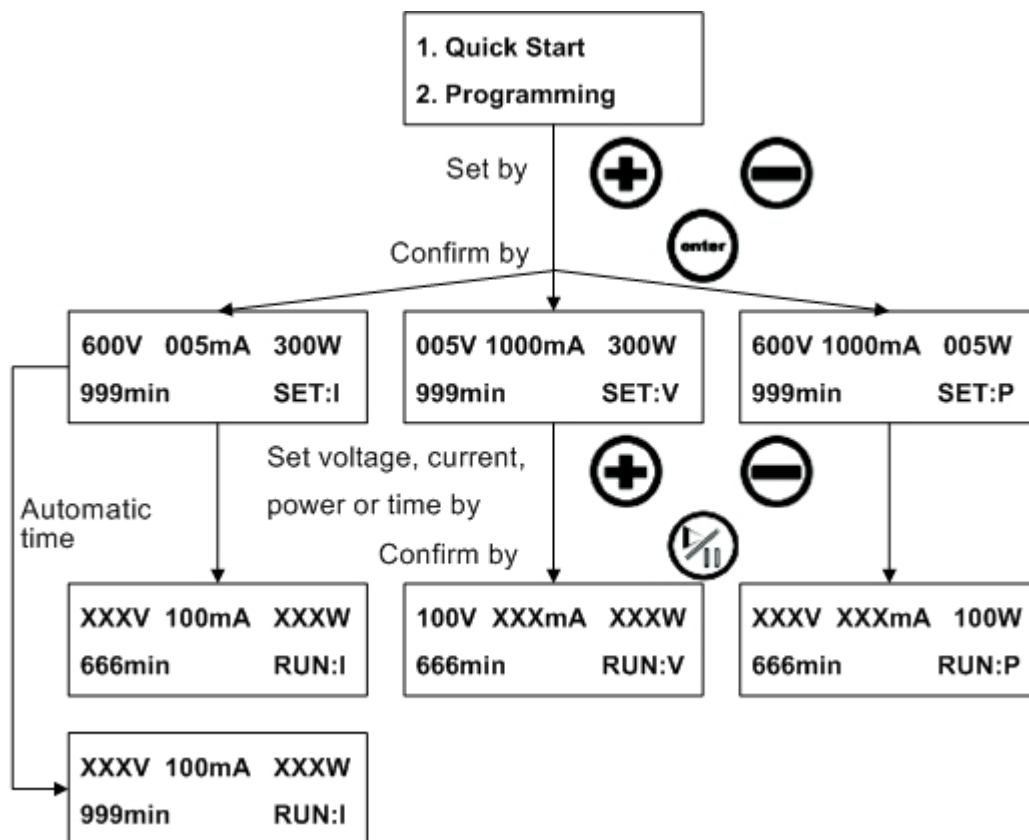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